

## *Sending your plots to the remote plotter*

### *Plotting on the large paper*

1. In the VNMR software, click on MAIN MENU, then on the second line at the far right (under the acquisition status window) click on MORE, then SELECT PLOTTER.
2. Each time you click SELECT PLOTTER, it will tell you the name of the selected plotter on the top of the page. It will continue to toggle through the list.
3. **plt1** - attached to the NMR for plotting      **prt1** – attached to the NMR for printing  
**remote-small** - the big printer on the table but 8½ x 11 paper  
**remote-big** – the big printer on the table but 11 x 17 paper
4. If you select anything but remote-big, the settings are done automatically.
5. If you select **remote-big**, you must make two changes...
6. On the command line type vp=12, this is a change in the vertical position of the spectrum
7. Click on MAIN MENU, then DISPLAY, then SIZE, then FULL. This will expand the spectrum to fit to the 11 x 17 size.